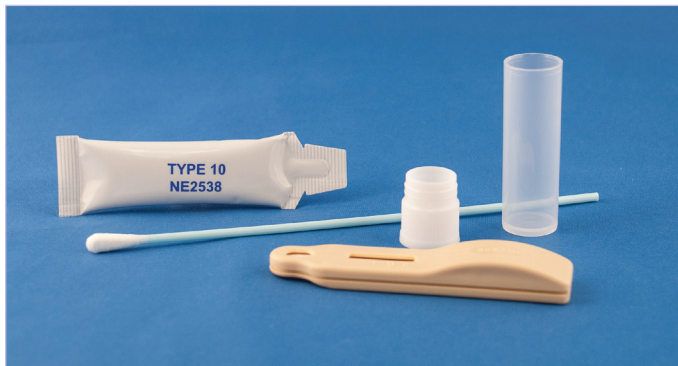


**Reveal® 3-D****For Sesame****Intended Use**

Reveal 3-D for Sesame is intended for the qualitative analysis of sesame residue in clean-in-place rinses and on environmental surfaces.

**The Test**

Reveal 3-D for Sesame is a unique hand-held device with a 3-line readout: a control line confirms the method has been performed successfully and two other lines differentiate nondetectable, low and high levels of sesame.

**The Procedure****Rinse and environmental samples:**

1. Add Type 10 buffer sachet to sample tube.
2. Add 1 mL (½ cap) of sample or environmental swab to a sample tube.
3. Cap and shake for 1 minute.
4. Remove cap and fill with liquid from tube.
5. Dip the Reveal 3-D device into the liquid in the lid.
6. Leave the cavity saturated until liquid is seen running into the test window.
7. Place device in a flat surface and allow test to develop for 5 minutes.

See package insert for complete instructions.

**Product Specifications**

Sensitivity:	Screens samples at 5 ppm sesame
Testing time:	5 minutes
Storage:	Refrigerated 2–8°C
Tests per kit:	10

**Results****Ordering Information**

Prod.#	Product description
8535	Reveal 3-D for Sesame